

B1 小動物室落菌數量

編號	區室	1月	2月	3月	4月	5月	6月	7月	8月	9月	10月	11月	12月
1	H20	1	0	0	0	2	1	1	0	0	0	0	0
2	H21	1	1	1	1	3	2	1	1	2	4	1	0
3	H22	1	1	1	1	4	1	3	6	0	1	0	0
4	H23	0	1	2	0	4	0	1	2	5	0	0	2
5	H24	0	2	0	1	6	1	2	3	3	3	1	3
6	H25	1	1	0	0	0	3	1	0	2	1	0	1
7	20-22 室走道	0	0	0	0	2	1	2	1	1	0	1	2
8	23-25 室走道	0	1	1	1	0	1	4	2	0	1	0	3
9	實驗室	1	1	0	1	5	2	1	1	1	0	0	0
10	洗衣室	0	1	0	2	0	1	1	5	3	1	1	4
11	滅菌室	0	0	1	0	0	1	0	1	0	0	1	0
12	清洗室	0	1	0	1	3	2	3	7	2	1	4	0
13	前走道	1	1	1	0	4	3	4	10	2	2	0	3
14	辦公室	1	1	1	0	1	3	3	6	2	9	4	0
15	更衣區	1	6	2	1	15	10	2	5	9	35	20	12
16	入口	0	3	1	1	6	2	4	1	11	7	6	2
17	20 操作台	0	0	0	0	0	0	0	0	0	0	0	0
18	21 操作台	0	0	0	0	0	0	0	0	0	0	0	0
19	22 操作台	0	0	0	0	0	0	0	0	0	0	0	0
20	23 操作台	0	0	0	0	0	0	0	0	0	0	0	0
21	24 操作台	0	0	1	0	0	0	0	0	0	0	0	0
22	25 操作台	0	0	0	0	0	0	0	0	0	0	0	0

1. BBL™ Stacker™ Plate TSB II， 5% Sheep blood (Becton, Dickinson)。
2. 方法：打開 Plate 上蓋，置於房間中央 30 分鐘，30 分鐘後蓋好上蓋，置入 37°C 培養箱中培養 48 小時，取出後計算落菌數量